Leica Cryofect

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Instructions for use

Keep only in original Container and up to 25 °C!

Highly flammable (VbF: B)!
Flash-point: 18 °C (DIN 51755)
Keep away from sources of ignition!
Do not spray into flames!
Attention with alcohol-sensitive surfaces!

Wear appropriate protective clothing when cleaning and disinfecting!
Disinfect the instrument and accessories every day!
Effectiveness of the disinfectant down to -20 °C experimentally tested.

1. Select a cryochamber temperature value down to -20 °C.
2. Remove the knife or blade from the knife holder.
3. Remove all samples, microscope slides and tools from the cryochamber.
4. Remove debris from the cryochamber.
Allow the cryochamber to reach the previously selected temperature.
Once the selected temperature is reached, either:
 - 5a. spray the disinfectant evenly on the contaminated surfaces - the surfaces should be covered with an even layer - or
 - 5b. soak a cloth with disinfectant and apply it on the contaminated surfaces.
6. Allow a reaction time of no less than 15 minutes.
7. Wipe it off with a tissue.
8. Dispose of tissue in compliance with the ruling waste disposal regulations of your institution.
9. Set the cryochamber temperature to the originally selected value.

If increased frost buildup occurs, start a manual defrost cycle.